gilt für / valid for
R 9000 + RS 9000

Bochem Instrumente GmbH
Laborbedarf

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www.bochem.de

Content

1 Introduction	12
2 Safety Instructions	12
2.1 Description of Symbols	12
2.2 Instructions	13
3 Product description	14
3.1 Key features	14
3.2 General survey	15
4 Packaging content	15
5 Positioning the device and commissioning	16
6 Job description	16
7 Maintenance and cleaning	17
8 Technical data	17
9 Certificate	18

1 Introduction

Thank you for purchasing our **Overhead Stirrer**. You did a good choice buying this product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Bochem reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of Symbols

WARNING

Information for avoiding injuries or fatal accidents.

ATTENTION

Safety instruction to avoid a damage of the product.

ATTENTION

Information about technical parameters.

Environment

Information about the operating environment.

WARNING of explosions

Safety instruction to avoid explosions.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- Do not use this device in areas with potentially explosive atmosphere.

- Make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- All circuits and electrotechnical components are **Bochem** owned technology. In fact of this it is necessary that only authorized personal repair the device. Contraventions exclude the warranty.

3 Product description


3.1 Key features

The **Overhead Stirrer** is designed for chemical, biological, medical and pharmaceutical applications.

Key features are as follows:

- Long life by simple and solid construction.
- Jog-Shuttle knob for easy and fast handling.
- Up to 3000 rpm.
- Max. viscosity: 10.000 mPas.
- Constant speed in spite of viscosity fluctuation.
- Motor provided with overload and overheat protection.
- Stepless variable adjustability of every stirrer concerning to the continuous bored chuck.
- Power bar graph and speed showed on illuminated LCD.
- Smooth increase of speed for higher safety.
- Storage function for speed value.

3.2 General survey

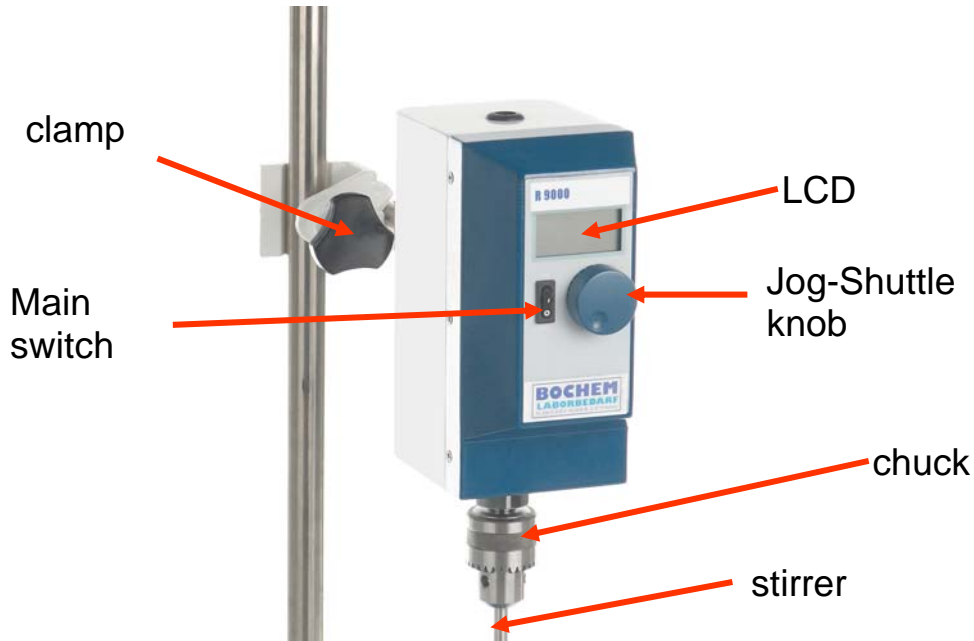4 Packing contents

quantity	R 9000
1	Allen key
1	body holder
1	chuck key
1	operating manual

quantity	RS 9000
1	impeller
1	stand & rod
1	Allen key
1	body holder
1	chuck key
1	operating manual
1	Clamp / Bosshead C 100

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Attach the clamp to the tripod rod.
5. Stick the device at the clamp and tighten the screw.
6. Check that each connection fits close.
7. Fix the stirrer in the chuck with included chuck key.
8. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
9. Make sure that you follow all safety instructions.
10. Connect the device with the power outlet.
11. Switch on the device and start the process.

6 Job description

- First switch on the device.
- If there is a value stored the stirring process starts by pushing the Jog-Shuttle.
- If there is no value stored the stirring process starts by turning the Jog-Shuttle.
- Change the set speed during the stirring application by turning the Jog-Shuttle.
- Start and stop by pushing the Jog-Shuttle.
- If you work with the same speed more often you can store this value and thereby you do not need to set often used speed values in every usage.
- To store values you only have to hold the Jog-Shuttle knob pushed for two seconds until the value in the first line in display is blinking.
- Delete stored values by over-storing with "0".

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull out the power plug of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you do not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Technical data

		R 9000
Stirring Capacity (H₂O)		Max.20 Lit.
Max. Viscosity (mPas)		10.000
Motor Rating (Input / Output)		130 / 120 W
Maximum Torque (Ncm)		145
Max. Shaft Diameter (mm)		10
Speed Range (rpm)		200 – 3.000
Controller		Digital Feedback Controller with Jog Shuttle knob (Turn + Push)
Display		Digital LCD Display (RPM and Power bar-graph)
Operating Environment	Ambient Temp. Range (°C)	0 ~ 40
	Max. Permissible Humidity (%)	80
Safety Device		Over Load & Heat Protector for motor
Dimension (wxdxh mm)		90x145x180
Weight (kg)		5,3
Others		Smooth Start and Change of the Stirring Speed Storage Function
Power Supply		220 V, 60 Hz

9 Certificate

Certificate

The quality and all features were checked by the manufacturer before the shipment.

We grant from the date of purchase

1 year guarantee

This certificate excludes damages caused by natural disasters or incorrect usage by the costumer.

Please complete the following table with the help of your account:

Artikel	Overhead stirrer R 9000
Typ	
Serial no	
Date	

Bochem Instrumente GmbH

Industriestrasse 11

D-35781 Weilburg

Tel: +49 (0)6471 / 92 82-0

Fax: +49 (0)6471 / 92 82-30

Email: info@bochem.de

www.bochem.de